

Liqui-Nox[®] Critical Cleaning Liquid Detergent



Use Instructions

Make a fresh 1% solution (2 1/2 Tbsp. per gal., 1 1/4 oz. per gal. or 10 ml per liter) in cold, warm or hot water. If available, use warm water. Use cold water for blood stains. For difficult soils, raise water temperature and use more detergent. Clean by soak, circulate, wipe or ultrasonic method. Not for spray machines, will foam. RINSE THOROUGHLY--preferably with running water. For critical cleaning, do final or all rinsing in distilled, deionized or purified water. For food contact surfaces, rinse with potable water. Used on a wide range of glass, ceramic, plastic and metal surfaces. Corrosion testing may be advisable.