

Luminox[®] Low-Foaming Neutral Cleaner



Use Instructions

For manual, soak and ultrasonic use, make a 3-5% solution (4-6 oz per gal. or 30-50 ml per liter) in warm or hot water. Protective gloves and eyewear are recommended. For spray machines, add detergent at the hot water wash cycle to make a 2% solution (2½ oz. per gal. or 20 ml per liter). For difficult soils raise water temperature and use more detergent. RINSE THOROUGHLY--preferably with running water. For critical cleaning do final or all rinsing in distilled, deionized or purified water. Corrosion testing may be advisable. Only the end user can make the determination as to whether this is appropriate or not.