

Diamond Knife Cleaning Instruction

Use Instructions



Fill the supplied beaker to 50 ml with a solution of 10% liquid detergent and water or other cleaning solution. Be sure to use only mild liquid detergent that do not contain ammonia or strong caustic chemicals, both of which can degrade the boat finish. Other recommended cleaning liquid are:

- A) 25% isopropyl alcohol in water or
- B) 25% acetone in water.

Clamp the diamond knife firmly in the holder and place it in the beaker. Set the timer to the desired time. Turn the cleaner "on" only after everything is in set up and ready to run. It will turn "off" automatically at the completion of the cycle (or manually by pressing the "off" button). The cycle time will vary from the indicated time if operating at 50 Hz.