


ZEROSTAT USE INSTRUCTIONS



SPI Supplies
206 Garfield Avenue,
West Chester, PA 19380, USA

Zerostat Antistatic Device #07008-AB

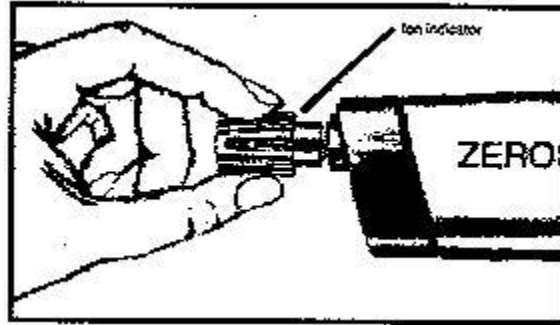

Hold the [Zerostat](#) within 12" (30.5 cm) of the surface or object to be treated. Squeeze the trigger slowly for about two seconds, to emit a stream of positive ionized air over the surface of the object. As the trigger is slowly released, a negative show of air ions is produced resulting in static neutralization. Repeat this procedure at several points across the surface of the object.

The Zerostat needs no batteries or electrical supply and unlike other antistatic products, the Zerostat contains no radioactive components whatsoever.

Important notice! Remove the ion indicator from the barrel before using the [Zerostat Antistatic Gun](#) to neutralize static charge. Ion indicator is for testing use only.

Caution! Do not use in an explosive or combustible environment.

11/17 EER